



## v-kit Pro Protocol Duplicator Modification

### **Background**


Up to and including v-kit Pro v 1.1.1 it has not been possible to duplicate previously duplicated protocols. This meant that simple changes to a duplicated protocol involved going all the way back to an original Supplied protocol and building a new protocol.

A new modification has been created (22-Aug-2011) which allows the duplication and publishing of protocols other than the original Supplied protocols.

### **How to use the protocol duplicator**

V:\Development\nlg10023 Protocol Duplication\vkitDupProtocolMods

Download then Unzip the folder named "vkitDupProtocolMods" This folder contains 11 files:

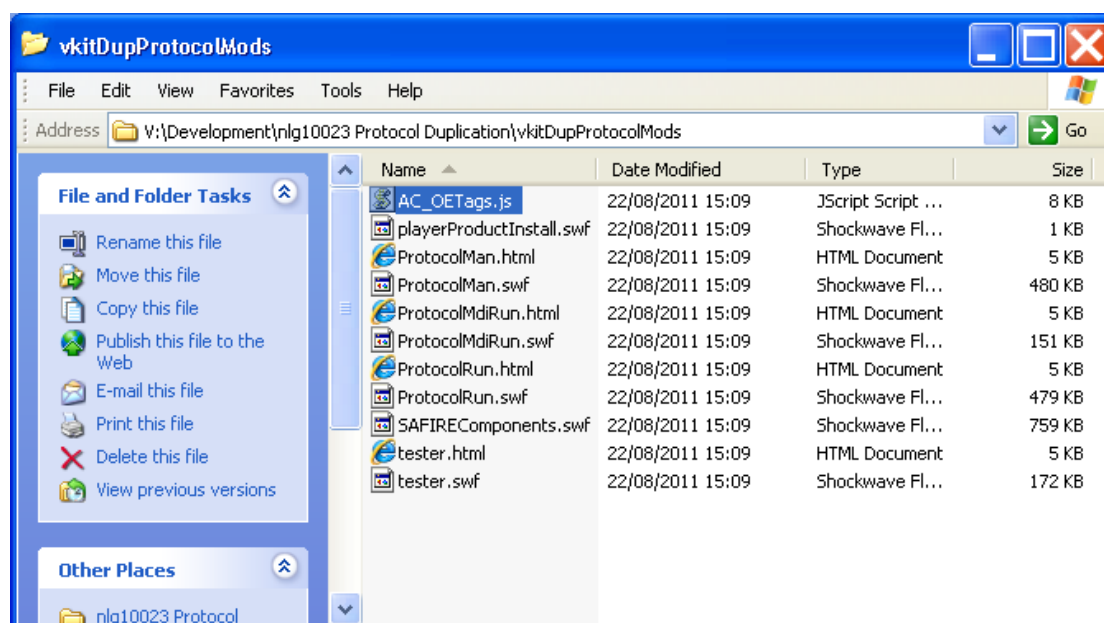

Locate your installation directory for v-kit Pro and drill down to the sub-folder named "docbase" i.e. C:\Program Files\v-kit Pro\dist\vkitbase\docbase

Make a copy your existing docbase folder and save somewhere safe as "docbase old" for example.

Copy the files you extracted previously into your docbase folder - it will overwrite all the older versions of the files (11 in total).

Run the v-kit application as normal. Verify that you can duplicate a non-supplied protocol, edit test defaults, re-name tests, publish the protocol and finally Run the protocol.

Neil Gowans, 09-Nov-2011